

Microcystins Strip Test Recreational Water 520022/520023

1. Collect Sample




Collect 1 to 2 mL of sample.


2. Transfer/QuikLyse™ *

Using the graduated pipette provided, transfer 1 mL of SAMPLE to the lysis vial containing the dried lysis reagent.

*QuikLyse™ reagents may be used in a method of U.S. Patent 9,739,777

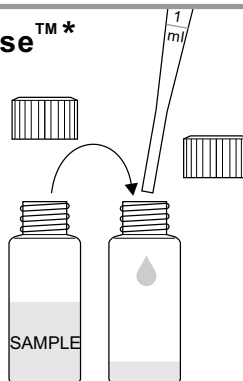

2 min.

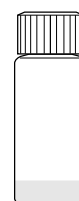



8 min.

Cap and shake for 2 minutes.

Let rest for 8 minutes.

3. Add Reagent Paper/QuikLyse™ *

Using the forceps provided, add 1 reagent paper to the lysis vial.

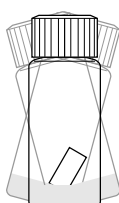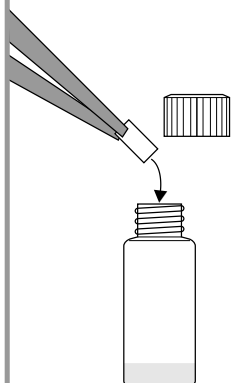

2 min.




8 min.

Cap and shake for 2 minutes.

Let rest for 8 minutes.

4. Transfer

Squeeze the upper bulb to draw sample into the pipette. The required amount of sample will fill the stem of the pipette. Any excess will flow into the overflow bulb. Squeeze the upper bulb again to dispense the required amount of sample into the conical tube.

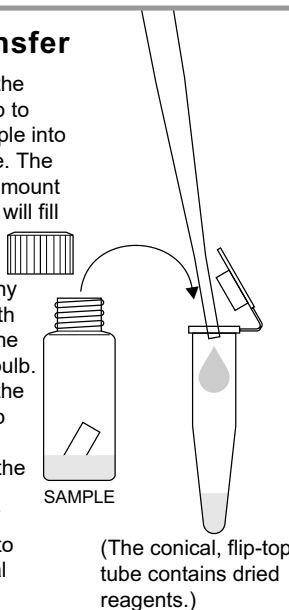

(The conical, flip-top tube contains dried reagents.)

5. Shake and Incubate



Close the conical, flip-top tube and shake for 30 seconds.



(Dried reagents will dissolve, turning the sample purple.)

6. Test

Insert test strip into conical, flip-top tube with arrow pointing down. (sample pad down).




Incubate for 10 minutes.




7. Dry

Remove test strip. Lay flat and allow to continue developing for 5 minutes.

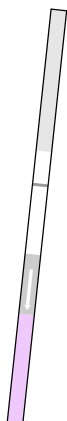

8. Interpret

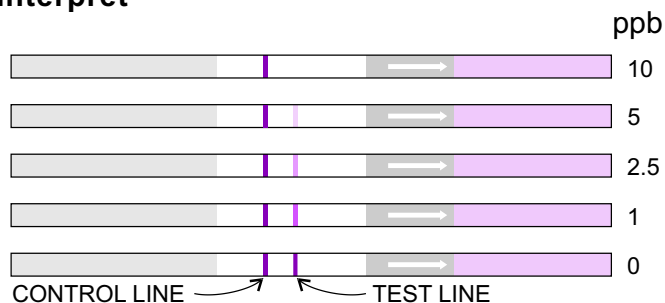

CONTROL LINE TEST LINE

INTERPRET TEST

CONTROL LINE	TEST LINE	INTERPRETATION
NO CONTROL LINE PRESENT	NO TEST LINE PRESENT	INVALID RESULT
CONTROL LINE PRESENT	NO TEST LINE PRESENT	>10 ppb
CONTROL LINE PRESENT	MODERATE INTENSITY TEST LINE PRESENT	BETWEEN 0 AND 10 ppb

Eurofins Abraxis
124 Railroad Drive
Warminster, PA 18974
WEB: www.abraxiskits.com

Date this Flow Chart is effective: 03FEB2022

T (215) 357 3911
F (215) 357 5232

Ordering: info.ET.Warminster@eurofinsus.com
Technical Support: support.ET.Warminster@eurofinsus.com

Version: 01